

**STANDARD OPERATING PROCEDURES**  
**DIVISION OF COMPARATIVE MEDICINE**  
**UNIVERSITY OF SOUTH FLORIDA**

SOP#: 1117.3

Date Issued: 11/00

Date Revised: 7/20

Page 1 of 2

---

|                        |                                                                                                  |
|------------------------|--------------------------------------------------------------------------------------------------|
| <b>TITLE:</b>          | <b>Refrigerator/Freezer Thermometer Use</b>                                                      |
| <b>SCOPE:</b>          | Research and Animal Care Personnel                                                               |
| <b>RESPONSIBILITY:</b> | Facility Manager, All Animal Program Personnel                                                   |
| <b>PURPOSE:</b>        | To Outline the Proper Procedures for Use and Maintenance of the Refrigerator/Freezer Thermometer |

---

**I. PURPOSE**

1. This procedure outlines the use and maintenance of a refrigerator/freezer thermometer used to monitor refrigerator/freezer temperature within the animal facilities.

**II. RESPONSIBILITY**

1. It is the responsibility of the Facility Manager to ensure that equipment is appropriately cleaned, maintained in good working order, and available for research personnel as requested.
2. It is the responsibility of the veterinary professional, administrative, and managerial staff to ensure that all research and technical staff using this equipment are adequately trained and experienced to perform hygrometer/thermometer determinations.

**III. EQUIPMENT USE**

1. Before using, slide open battery cover located on back of unit and insert one AA battery. Replace cover.
2. Slide the switch on back of unit to °C for Celsius.
3. Slide alarm to off. Refer to instruction booklet if necessary to use the alarm mode.
4. Insert external sensor probe jack into outlet on side of unit.

8. Press Memory clear button to reset the MAX/MIN temperatures.

#### IV. MAINTENANCE

1. Clean unit by wiping with a mild disinfectant and soft cloth.
2. Replace AA alkaline battery when digits flash, there is no display, or display is faint.
3. Abnormal readings not corrected by battery replacement should be promptly reported to the Facility Manager.
4. Annually recertify and label the unit in accordance with **SOP #1115**  
***Thermometer/Hygrometer Certification.***
4. Any additional maintenance/service should be performed by authorized personnel.

#### V. REFERENCES

Manufacturer's Manual.

Approved:

Date: